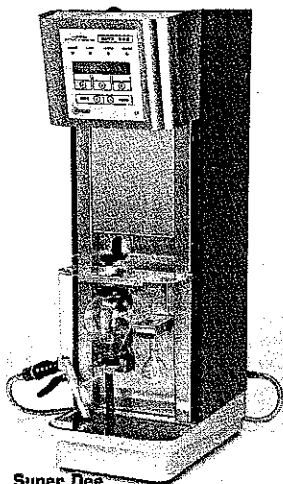
Misure automatiche a norma

Super Dee

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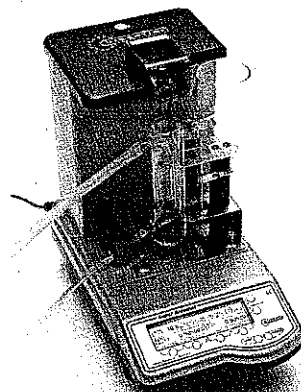
Super Alcoma

Automatic measurements complying with law

An oenological distilling unit and an electronic hydrostatic balance enable to make fast and accurate analysis, complying with the community rules in force

Gibertini's new digital distilling unit Super Dee complies with the official method for the determination of the alcoholic strength by volume in wines, musts and alcoholic. Its execution times for 100 ml wines accounts for 3-4 minutes, for 200 ml wines for 6-7 minutes; volatile acidity, 250 ml 6-7 minutes, while for alcoholic beverages: 4-8 minutes. Control panel features easy-to-read alphanumeric intelligent display and keyboard. It drives and controls the microprocessor for: the regulation and the tare of the new balance, the volume to be distilled, the real time value of the distilled amount; check of the inlet cooling water: temperature, pressure and flow, the different languages, the warning of missing glassware, the warning of opened protection or drainage. No more distilled water is required due to the integrated steam generator. Direct connection to the tap water and cartridge for demineralisation interchangeable by the user. Available also without steam generator. The following new technical characteristics have been introduced: new vertical sliding plexiglass protection; new phial with enlarged drainage to distillate wine lees and grape pulp; new system rapid and operational (one hand operation) for

holding the glassware. It inhibits the distillation if the glassware is not present; new shower for washing the phial with flow regulation; magnetic filter at the inlet of the water for stopping heavy particles; automatic stop of the instrument after the distillation, no wasting of water. The Superalcomat electronic hydrostatic balance stands out for the automatic determination of the TAV, the total dry extract and musts evaluation. The readout of the alcoholic strength by volume is shown on a large LCD screen simultaneously with the readout of temperature, density at 20 °C and other parameters. The reading module is provided with a shower proof keyboards whose functional keys allow: correction of the floater volume with H2O or certified standard solutions, zeroset and calibration of the super Alcomat, date, hour, minute setting and other functions; the printing of the reading values and other parameters (printer not included), must evaluation, Baumé, Babo, Brix and Dechsle degrees; working temperature: 10-30 °C optimum 15-25 °C, languages. The special microprocessor drives the functions and makes the internal problem diagnostics for preventive maintenance.