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*All the information given in this manual is what available at the time of printing.
The manufacturer reserves the right to make changes to the product at any time without notice.
It is advisable to verify if there are any updates.*

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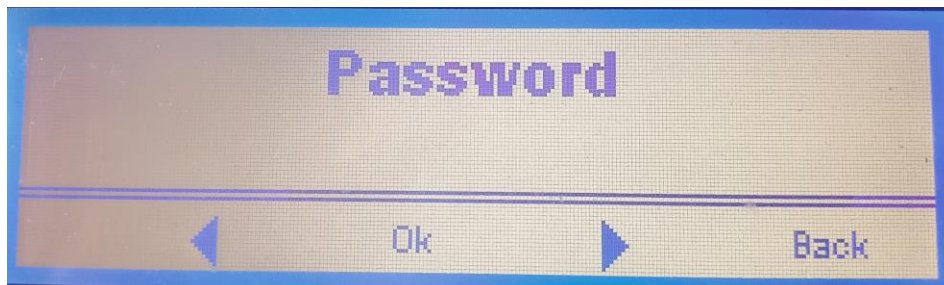
All rights reserved.

*Forbidden **any form** of print, duplication, reproduction and publication
of this manual, or part of it, without the written agreement of
GIBERTINI ELETTRONICA Srl*

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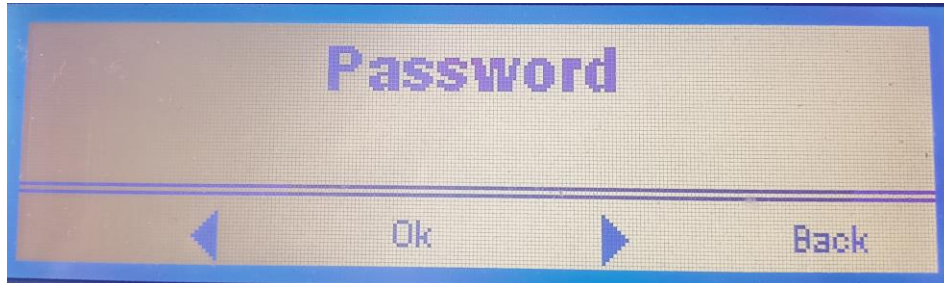
Password management

From the Setup Parameters menu select Password and confirm with "ok"

With the + and - keys change the selected value and with next go to the next digit. It is possible to confirm with "ok" or "Back" to move backwards or cancel the operation with the "tare" key.

Once the password has been enabled, it will be necessary to enter it in order to access:

- Setup parameters
- Date & Time Adjust
- External calibration

The new scale has a default password of '0000' which corresponds to a disabled password, once a password has been entered.

To disable the password protection: log in with the current password and type '0000'

If you forgot the current password : it is possible to reset the password by restoring the factory data, this operation will not return the value to '0000' but will use the first four digits of the instrument's serial number, for example if the serial number is ' 156787 'after a factory reset the password will still be enabled and will have the value ' 1567 '

info@gibertini.com
Tel. (+39) 023541434

ISO 9001:2015

LAT N° 094

GIBERTINI
ELETRONICA s.r.l.
Via Bellini 37
20026 Novate Mil.se
Milano (Italy)

Tel. (+39) 02 3541434
Fax (+39) 02 3541438
www.gibertini.com
sales@gibertini.com