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*All the information given in this manual is what available at the time of printing.  
The manufacturer reserves the right to make changes to the product at any time without notice.  
It is advisable to verify if there are any updates.*

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*All rights reserved.*

*Forbidden **any form** of print, duplication, reproduction and publication  
of this manual, or part of it, without the written agreement of  
GIBERTINI ELETTRONICA Srl*

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**NOTE**

Read the instructions carefully before operating the Balance.  
This is an accuracy instrument and must be handled with care.

**INDEX**

<b>1 WARNING</b>	<b>23</b>
<b>2 TECHNICAL CHARACTERISTICS</b>	<b>24</b>
<b>3 LOCATION</b>	<b>24</b>
<b>4 INSTALLATION</b>	<b>25</b>
<b>5 OPERATING CONTROLS OF THE BALANCE</b>	<b>26</b>
<b>6 OPERATING INSTRUCTIONS</b>	<b>26</b>
<b>7 KEYBOARD FUNCTION</b>	<b>26</b>
<b>8 TARE</b>	<b>27</b>
<b>9 MAIN MENU AND OPERATING FUNCTIONS</b>	<b>27</b>
9.1 Instrument Functions	28
9.1.1 Simple Weighing	28
9.1.2 Piece Counting	28
9.1.3 Percentage	29
9.1.4 Standard Deviation	29
9.1.5 Density	29
9.1.6 Dynamic Animal Weighing (DYNAMIC WEIGHING FUNCTION FOR ANIMALS)	30
9.2 2nd Unit	31
9.3 Setup parameters	32
9.3.1 Serial Port	32
9.3.1.1 Baud Rate	32
9.3.1.2 Byte Format	33
9.3.1.3 Protocol	33
9.3.1.4 Transmit Option	33
9.3.1.5 Stable Option	33
9.3.2 Printer Type	33
9.3.3 Calibration	33
9.3.4 Auto Zero	34
9.3.5 Filtering	34
9.3.6 Idle Setting	34
9.3.7 Lcd Contrast	34
9.3.8 Time & Date Format	34
9.3.9 Languages	34
9.3.10 Password	34
9.3.11 Set Default	35
9.4 Date & Time Adjust	35
9.4.1 Time Setting	35
9.4.2 Date Setting	35
9.4.3 Day of Week Setting	35
9.5 WiFi	35
9.6 External Calibration	37
<b>10 QUICK RESTORE OF DEFAULT PARAMETERS</b>	<b>37</b>
<b>11 SERIAL OUTPUT</b>	<b>38</b>
11.1 SERIAL OUTPUT CONNECTION	38
11.2 DATA TRANSMISSION FORMAT	38
11.2.1 CRYSTAL protocol	38
11.2.2 EURO protocol	39
11.2.3 PRINTER protocol	39
11.3 SERIAL COMMANDS	39

## 1 WARNING

The power supply pack of the EU-C Gold scales complies with Electrical Protection Class II. Before connecting the power, check that the operating voltage, indicated on the Power supply pack, matches the line voltage (100 / 240 Vac).

**EU-C GOLD SCALES MAY ONLY BE USED WITH THE POWER SUPPLY PACK.**

**THE BALANCE MUST BE OPERATED ONLY IN DRY, CLEAN AND IF POSSIBLE VIBRATION-FREE ENVIRONMENT.**

**THE BALANCE MUST NOT BE OPERATED IN AREAS WHERE THERE ARE EXPLOSION RISKS.**

### NOTE:

In case of improper setting of the balance parameters during setup, the balance may power-up in an incorrect configuration.

The user can restore the Factory configuration to avoid the malfunction as follows:

Disconnect the balance supply.  
Keep the first key on the left pressed  
Reconnect the supply  
Wait for display showing 'Set Factory Parameter'  
Release the key  
Now the Factory default settings are restored.

**ALL ELECTRONIC BALANCES, WHICH ARE SENSITIVE TO CHANGES OF GRAVITY, MUST BE CHECKED AND CALIBRATED IN THE PLACE WHERE THEY ARE TO BE USED (CE 2014/30/UE).**

**ALL OUR INSTRUMENTS MEET INTERNATIONAL EMC-EMI-RFI STANDARDS ACCORDING TO CE 2014/31/UE DIRECTIVE ON ELECTROMAGNETIC COMPATIBILITY.**

## 2 TECHNICAL CHARACTERISTICS

Model - EU-C GOLD	802	2002	4002	4000	7500PQ	7500PT	7500DR	10000PT
Capacity	820g	2000g	4200g	4200g	7500g		950 / 7500	10200
Readability	0.01g	0,01g	0,01g	0,1g	0,1g		0.01 / 0.1	0.1
Linearity	± 0,05	± 0,02	± 0,02	± 0,3	± 0,3		± 0,05 / ± 0,3	± 0,5
Repeatability	± 0,02	± 0,01	± 0,02	± 0,2	± 0,2		± 0,02 / ± 0,2	± 0,2
Response time (seconds)	2	2	2	1	1		2 / 1	2
Pan dimension (mm)	150x150	150x150	150x150	Ø 190	190x190	Ø 190	Ø 190	Ø 190
Warm-up time (minutes)	≥ 15							
External calibration (ClasseF1) multiples of	100g	1000g	1000g	1000g	1000g	1000g	1000g	1000g
Operating temperature	10 ÷ 40°C							
Data output	RS232 I/O - USB 2.0 - WiFi							
Operative voltage	100 / 240 V ac -15/+10% (from 85 to 264 V) 50/60Hz							
Power consumption mA	200							
Dimension	216 x 320 x 100 mm (W x D x H)							
Weight	3 Kg (power supply included)							
Standard equipment	Power supply - User manual - CD with driver							

## 3 LOCATION

Choose an appropriate location taking into account these criteria:

- ◆ Vibrations and air drafts in the ambient must be avoided.
- ◆ The supporting stand must be rigid.
- ◆ The main line must be stable. It is not advisable to use the same electric line of the fluorescent lamps or other electric apparatus that can be switched on and off randomly.
- ◆ The balance must be kept far from magnetic fields.
- ◆ Environmental conditions with wide temperature changes must be avoided. Ambient temperature above 30 °C or below 10 °C may affect the operation of the balance and its accuracy.

In very dry environments electrostatic charges may build up, especially if glass or porcelain objects are used


## 4 INSTALLATION

### Instructions to remove the locking device of the balance

Unpack the balance and all its accessories.

- ◆ Unscrew the knob on the locking bracket as indicated by the arrow (anti clockwise).  
Retain the locking bracket in case the balance has to be returned to the supplier. This bracket can be refastened to the balance by replacing the knob.
- ◆ Screw the pan with the central screw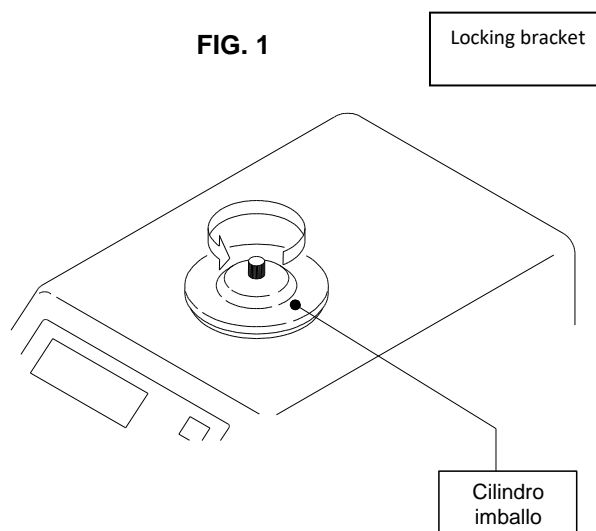

- ◆ Level the balance using the spirit level on the rear and rotating the levelling feet under the front part of the basement (Fig 2).
- ◆ Insert the power supply cable located on the rear panel (Fig 2).
- ◆ Insert the power supply in the main line plug checking the supply voltage printed on the power supply case (100 / 240 Vac ) - (Fig 2).

## 5 OPERATING CONTROLS OF THE BALANCE

1. Multifunction keys
2. Tare Key
3. Graphic Display

## 6 OPERATING INSTRUCTIONS

Wait for about 60 minutes after power up before performing any weighing operation. It's advisable to leave the balance permanently connected to the Ac line.

## 7 KEYBOARD FUNCTION

EU-C Gold series scale keyboard has a 4 multifunction keys and a Tare key. In some cases, depending on the activated menu, multifunction keys, if pressed for less than two seconds, activate the first function, if pressed for more than two seconds they activate the second function, as indicated in the respective specific paragraphs.

## 8 TARE

After a container is placed on the pan, press the “T” (TARE) key.

The LCD display will show zero value.

The balance has a tare capacity up to its measuring range.

Place the sample to be weighed at the center of weighing pan.

The display shows the sample net weight.

It's possible to use the tare function after adding several samples up to the maximum balance capacity indicated by a bargraph (in graphic mode and % of full scale) on the left side of the LCD display.

## 9 MAIN MENU AND OPERATING FUNCTIONS

The balance can be operated in the following configurations:

Balance with gram, carat, ounce, pound measuring unit

Piece counting balance

Percentage computing balance

The selection is performed as follows:


Hold MODE key pressed until the display shows “UNIT 2”, release the key. Pressing key MODE (short time) it is possible to select the functions “UNIT 2”, “SETUP”, “CAL E”, wrapping around at the end of the list.

The ON/OFF key activates the selected submenu, while the TARE key exits the main menu.

When the submenu is entered, the keys operate as above: pressing ON/OFF the selection is confirmed, while pressing TARE the selection is left unchanged.

NOTE: If no key is activated for 20 seconds, the menu level is incremented up to the end, without altering any parameter and the balance resumes its normal operation.


**Fig. 5 – Graphic display detail**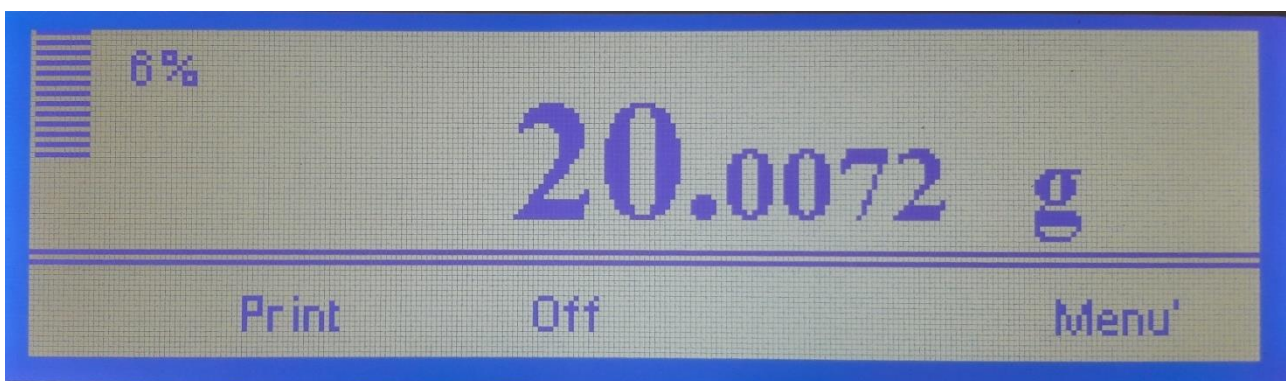

- When entering a menu or sub-menu, the Back key allows you to return to the previous menu
- Within each menu, if no button is pressed for about 20 seconds, the main menu is displayed again.
- Within the menus subject to modification, any modified parameter always requires confirmation with the OK button

### Main Menu Flow


To enter the main menu, press the Menu button for at least 2 s  
 The following continuous scrolling submenus are displayed: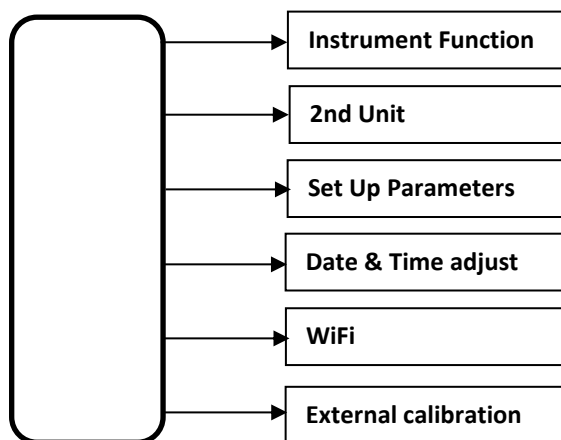

To enter the submenus, scroll with the arrows until the desired submenu is displayed and confirm with the OK button

## 9.1 Instrument Functions

To enter the Instrument Function menu, press the Menu button for at least 2 s and confirm with OK  
 The following continuous scrolling submenus are displayed: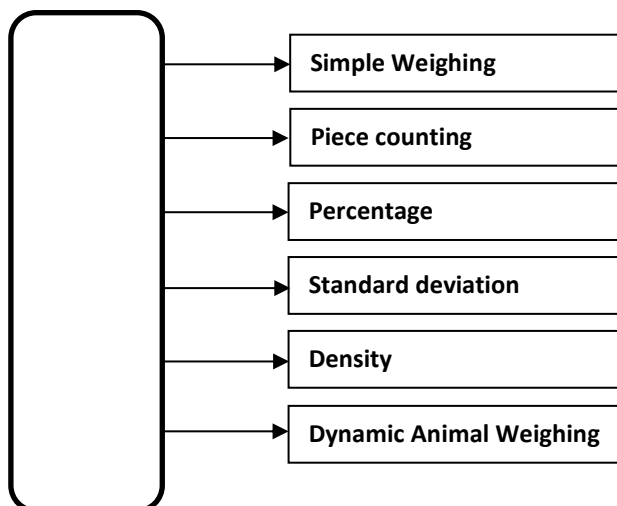

### 9.1.1 Simple Weighing

Press the Simple Weighing button to exit the Instrument Functions submenu or to disable any active Instrument Function

### 9.1.2 Piece Counting

To activate the Piece Counting function, press the OK button from its submenu.

It is displayed the screen from which it is possible to select, by pressing the Sample button, the quantity of

pieces to be weighed by sampling among the values 5, 10, 20, 50, 100.

When the number of pieces to be sampled has been placed on the pan and the stable data symbol is lit, press the Acquire key to confirm.

The scale automatically calculates the unit weight and starts weighing with the piece-counting function.

**NOTE:** In case of sampling with too low values, the system signals an error and the display shows the wording PCS ERROR.

To exit the error condition, press the T key.

### 9.1.3 Percentage

To activate the Percentage function, press the OK button from its submenu.

It is displayed the screen from which it is possible to select, by pressing the Percentage button, the storage mode of the 100% value.

Place the reference sample on the weighing pan.

When the stable data symbol is lit, press the Percentage key to confirm by weight corresponding to 100%.

The scale acquires the reference value and starts weighing with a percentage calculation function.

All the weighs that will be carried out after the acquisition will have the sample previously acquired as a percentage reference.

They appear in two sub-fields:

- 'T' Target the value corresponding to 100%.
- 'S' Sample the value of the sample under examination.

There are two modes of indication of the percentage, using the Mode n button it is possible to switch between the two operating modes. In mode 1 the percentage considers the target as 100%, in mode 2 the target is considered 0% and the percentage variation + or – with respect to the target is indicated.

EU-C Gold can measure percentage values of 500% with respect to the sample, provided that the maximum capacity of the scale is not exceeded.

**NOTE:** In case of sampling with too low values, an acquisition error will be signaled and the display will show the message PERC ERROR.

To exit the error condition, press the T key.

### 9.1.4 Standard Deviation

To activate the Standard Deviation function, press the OK button from the relevant submenu.

It is displayed the screen from which it is possible to select, by pressing the Deviation key, the method for storing the weighing values that will determine the calculation of the standard deviation.

To store the chosen weight value to carry out the weighing in sequence, press the Acquire key once the stable weight condition has been reached.

Each weigh made is stored with a progressive number of "Samples".

The calculation of the Standard Deviation is performed and shown on the display in real time starting from Sample n. 2.

In the event of a weighing error, it is possible to delete the last data acquired by pressing the Clear Last button.

At the end of the acquisitions, pressing the Results button displays the result screens which include the following data: number of weighs (Sample), average value of weighs (Average), Standard Deviation (Deviation), Total sum of weighs (Total Sum), weighing with minimum value (Lowest), weighing with maximum value (Highest), Difference between weighing with minimum and maximum value (Max-Min). By pressing the Print button it is possible to print the receipt which includes all the data listed above.

### 9.1.5 Density

To activate the Density function, press the OK button from the relevant submenu.

It is displayed the screen from which it is possible to select, by pressing the Density key, the method for storing the weighing values that will determine the density calculation.

The interactive menu allows to guide the operator in storing the values necessary for the calculation through the following messages on the display:

- WEIGHT IN AIR: Place the object to be weighed on the pan, wait for the stable weight condition and store the value by pressing the Acquire key
- WEIGHT IN LIQUID : place the object to be weighed on the weighing device in the liquid, wait for the stable weight condition and store the value by pressing the Acquire key
- The display shows the density value relating to the two stored weighs

The density calculation is performed considering the density of water at 20°C (0.998205) if you use a different liquid and/or work at a different temperature by pressing the Liquid dens key for at least 2 seconds you can access the input function of an arbitrary liquid density, **N.B. it must be the user's responsibility to check that this value corresponds to the density of the liquid at the current working temperature.**

**To proceed with a new density determination, press the Exit button and then Density.**

### **9.1.6 Dynamic Animal Weighing (DYNAMIC WEIGHING FUNCTION FOR ANIMALS)**

To activate the Dynamic Animal Weighing function, press the OK key from the relative submenu.

The screen is displayed from which it is possible to select, by pressing the Dynamic key, the sampling time of the weigh among the values of 1, 2, 4, 6.8, 10, 15, 20 seconds.

When the acquisition time has been chosen and the animal to be weighed has been placed on the pan, press the Acquire key to confirm.


The scale automatically starts acquiring the weighs displaying the residual time, after which the message ACQUIRED and the average value of the readings will be displayed.

TO EXIT THE "INSTRUMENT FUNCTIONS" SUBMENU PRESS THE "Simple Weighing" KEY

## 9.2 2nd Unit

To activate the 2nd Unit function, press the OK button from the relevant submenu.


The current selected unit of measurement is displayed and the alternative units can be selected using the two arrows. To store the desired unit of measurement, press the OK button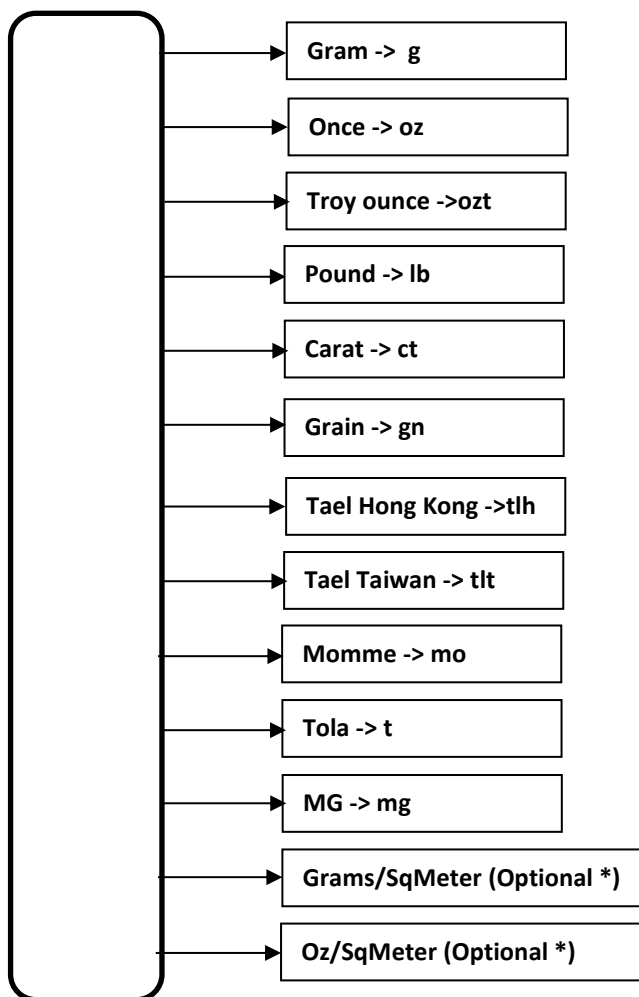

\* Optional units of measure that require the use of the appropriate sample cutting tool not supplied with the balance.


### 9.3 Setup parameters

To enter the Setup Parameters submenu, press the OK button from the relevant submenu  
The following continuous scrolling submenus are displayed: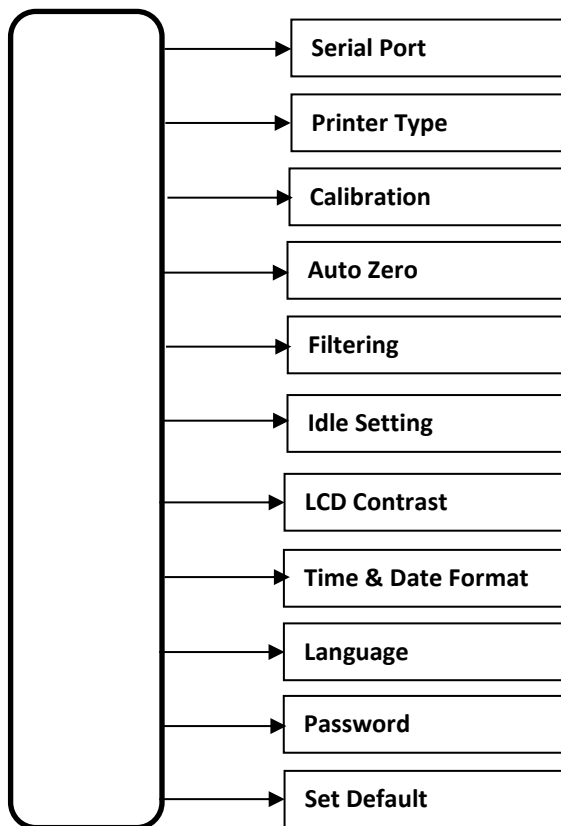

#### 9.3.1 Serial Port

To enter the Serial Port submenu, press the OK button from the relevant submenu  
The following continuous scrolling submenus are displayed:

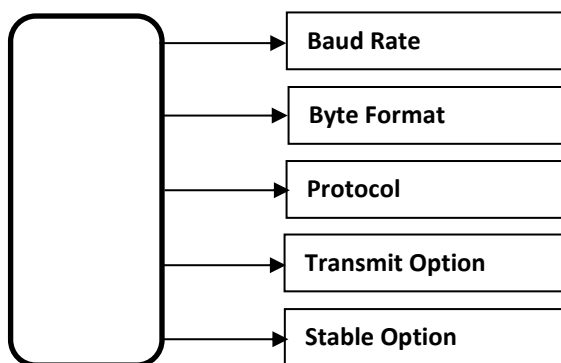##### 9.3.1.1 Baud Rate

To enter the Baud Rate function, press the OK button from the relevant submenu  
To change the serial transmission speed select between 1200, 2400, 4800, 9600, 19200 and confirm with the OK button

### 9.3.1.2 Byte Format

To modify the serial transmission data format select between 7E1, 7O1, 7N2, 7E2, 7O2, 8N1, 8E1, 8O1, 8N2 (as indicated in the following table) and confirm with the OK key

DATA FORMATS	DATA BIT	PARITY CHECK TYPE	BIT STOP
7 E 1	7	EVEN	1
7 O 1	7	ODD	1
7 N 2	7	NO PARITY	2
7 E 2	7	EVEN	2
7 O 2	7	ODD	2
8 N 1	8	NO PARITY	1
8 E 1	8	EVEN	1
8 O 1	8	ODD	1
8 N 2	8	NO PARITY	2

### 9.3.1.3 Protocol

To change the serial data output protocol select from Europe, Crystal, Printer and confirm with the OK button. Refer to the DATA TRANSMISSIONS section for more information.

### 9.3.1.4 Transmit Option

To change the serial data transmission mode, select between Request and Continue and confirm with the OK button.

**Request** The data is transmitted only if there has been a request sent to the serial port

**Continue** The data is transmitted continuously at each cycle, even if there has been no request from the serial port. It is useful when the scale must transmit data to a peripheral that is unable to send commands to the scale

### 9.3.1.5 Stable Option

To change the transmission mode to stable data or not, select between Stable Mode OFF and Stable Mode ON and confirm with OK. In ON mode, the data is transmitted to the request only if the weight value on the pan is stable, otherwise it waits for the stability of the data and then transmits. In OFF mode the data is transmitted independently of the weight value stability.

### 9.3.2 Printer Type

To change the type of printer connected to the serial, select from Thermal Printer (24 column), Thermal Printer (40 column), Impact Printer (24 column), Impact Printer (40 column) and confirm with the OK button.

### 9.3.3 Calibration

To modify the mode in which the calibration process is carried out select between Calibration AUTO, Calibration OFF, Calibration ON and confirm with the OK key.

The calibration mode can be selected by the user according to the following:

<b>Calibration OFF</b>	No calibration process can be performed. By selecting this option, the possibility of external calibration CAL E and internal (if present) are inhibited.
<b>Calibration ON</b>	With this option active, it is possible to carry out with the relative commands both external and internal calibration (if present). On the other hand, timed internal calibration is not enabled.

### 9.3.4 Auto Zero

To change the mode in which automatic zero tracking is performed, select between Zero ON, Zero OFF and confirm with the OK button.

**ZERO ON** If the Auto-Zero function is active and if there is no mass on the pan, the zero tracking algorithm keeps the value measured at zero. The function is useful for canceling small zero drifts.

**ZERO OFF** With the Auto-Zero function in the OFF position, zero correction is not performed.

### 9.3.5 Filtering

To change the dynamic behavior of the scale, select from Slow Filtering, Average Filtering, Fast Filtering (Dosage) and confirm with the OK button.

**Slow Filtering:** slow response time. Suitable for environments disturbed with air currents and / or vibrations

**Average Filtering:** average response time between slow and fast. Suitable for general work environments.

**Fast Filtering:** fast response time. Suitable for acquiring rapid weight changes (e.g. dosages).

### 9.3.6 Idle Setting


To change the mode in which the scale automatically switches off the display following non-use for a specified time, select between **Idle OFF** (no auto power off), **Idle ON** (auto time off) and confirm with the OK key.

### 9.3.7 Lcd Contrast

To adapt the brightness of the display to the work environment, select the function and use the  $\pm$  keys to adjust the intensity in% ( $\pm 5\%$ ), confirming the preferred selection with the Update key.

### 9.3.8 Time & Date Format

To change the way the balance displays Date and Time, select the options indicated in the following diagram and confirm with the OK key.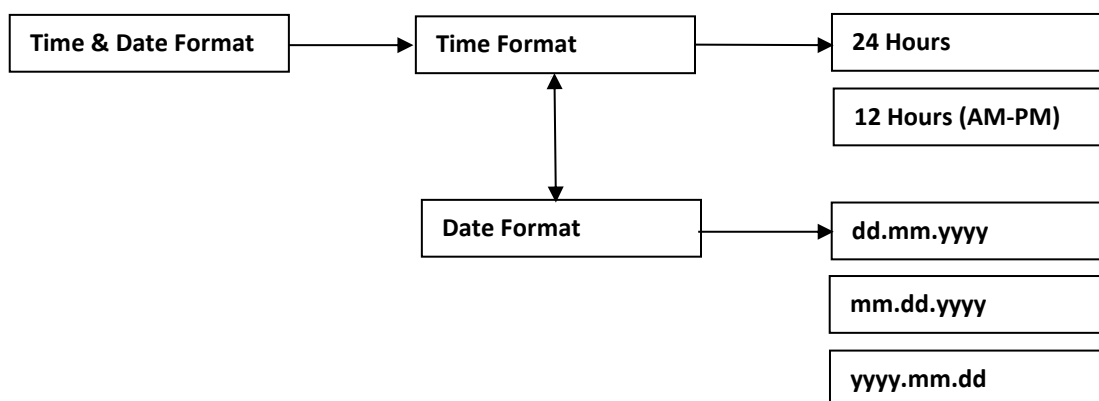

### 9.3.9 Languages

To change the display language, select from English, Italiano and confirm with the OK button.


### 9.3.10 Password

Enter a password to protect the functional parameters of the instrument.

You can use a 4-character numeric password, the default value is '0000' which corresponds to a disabled password.

From the Setup Parameters menu select Password and confirm with "ok"

With the + and - keys change the selected value and with next go to the next digit. It is possible to confirm with "ok" or "Back" to move backwards or cancel the operation with the "tare" key.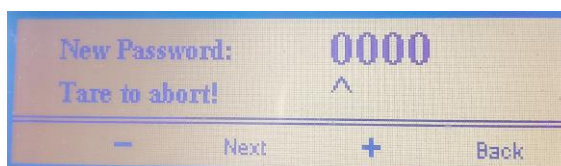

Once the password has been enabled, it will be necessary to enter it in order to access:

- Setup parameters
- Date & Time Adjust
- External calibration

For password management details see attachment.

### 9.3.11 Set Default

All the parameters modified in the SETUP menu are immediately operational as soon as they are changed. However, if the scale is turned off by removing the power supply, they will be lost.


As a consequence of this, the next time the power is turned on, the scale will work with the previous set parameters.

In order to maintain a personalized configuration, it is possible to store each change by selecting **Store parameters & exit** or **Exit without saving** option to not store the changes made. Confirm your selection with the OK button.

## 9.4 Date & Time Adjust

To enter the Date & Time Adjust submenu, press the OK button from the relevant submenu

The following continuous scrolling submenus are displayed: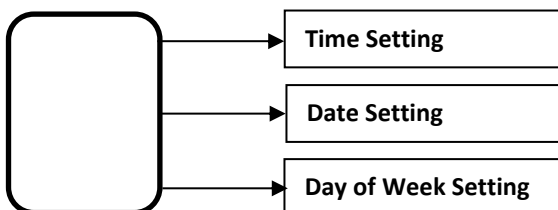

### 9.4.1 Time Setting

To adjust the hour displayed on the scale, enter the relative submenu with the OK button and adjust the hours and minutes using the  $\pm$  and Next buttons. Pressing Next and Update successively the new time is stored.

### 9.4.2 Date Setting

To adjust the date displayed on the scale, enter the relevant submenu with the OK button and adjust the day, month and year with the  $\pm$  and Next buttons. Pressing Next and Update successively the new date is stored.

### 9.4.3 Day of Week Setting

To adjust the day of the week displayed on the scale, enter the relative submenu with the OK key and use the  $\pm$  keys to select the day of the week, day, month and year. Pressing Update later the new day of the week is stored.


## 9.5 WiFi MENU

For the configuration of the network parameters it is not necessary to use any dedicated software, the whole procedure can be performed with a normal internet browser (Firefox, Chrome, Edge etc.) and from any device and / or operating system, simply connect to the configuration page with this procedure:

- Open the search for wireless networks, the factory configuration of the wireless network is unique for each instrument and is composed as follows:

- **SSID** = EU-C\_XXXXXXXXXX (where XXXXXXXXXXXX correspond to the MACADDRESS of the device)
- **PASS** = EU-C-PASS


- in the example image, the network has SSID EU-C\_3C71BF2F9CAC:

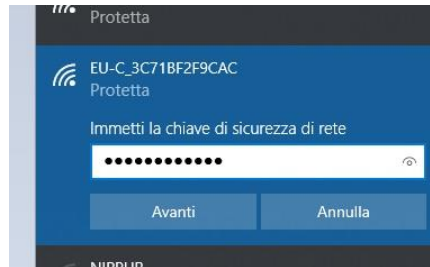
- Once connected to the network by accessing with a browser at 192.168.4.1, a summary page of the current configuration will be displayed:

- By accessing the setup page, it will be possible to configure the instrument in Access Point or Station mode.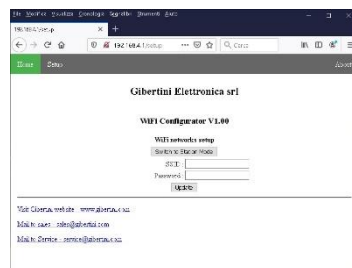

- In Access Point mode enter the desired name for the network and the relative password, in this mode the scale is independent and will create the required network autonomously without the need for any network access point. The network address in this mode will always be 192.168 .4.1, it will then be possible to use the WiFi connection with the scale even in the absence of an existing network.
- In station mode, the scale will scan the networks present in the work site, it will therefore be possible to select your work network from the list by entering the correct configuration parameters:
  - **IP Addr:** Enter a valid IP address for your network or leave DHCP to get one automatically.
  - **Password:** Enter your network password if present or leave blank for open network. (n.b. the networks detected will be displayed in the list with the final '\*' if password protected).
- Once the configuration has been complete, the scale will start the connection procedure to the local network, showing "Connecting" on the display. If the connection is established, will show "Connected" on the display for about 1 second and the new IP address will follow. The brief summary page of the new configuration will be displayed in the browser.

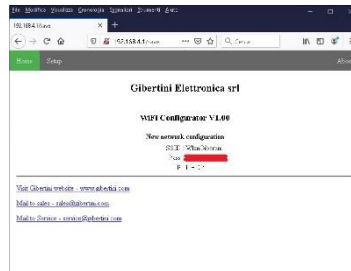- If the connection is refused, the display will show "Error" for about 2 sec. The access point functionality will be restored and will be displayed the default IP address 192.168.4.1

Once the configuration is complete it will be possible to communicate with the scale via TCP protocol on port 333, according to the standard communication protocol normally used for the RS232 connection.

## 9.6 External Calibration

Scales with a capacity of 4200 g, 6300 g and 10200 g use a mass of 1000 g as the first calibration weight; scales with a capacity of 2200 g use a 500 g mass as the first calibration weight; the 210 g scale uses a 100 g mass as the first calibration weight.

By activating the External Calibration function, a warning message indicates the indications relating to the masses to be used for external calibration. See also Chapter 2 - Technical characteristics

To enter the External Calibration submenu, press the OK button from the relevant submenu

When this function is activated, the calibration process is immediately performed.

Automatically you exit all menu functions and the message "Phase 1 Load weight" appears on the display.

In this phase, the scale waits on the weighing pan for a calibration mass whose weight must be a function of the full scale value of the scale itself.

EU-C Gold automatically checks and verifies the calibration weight and accepts multiple values of the base weight of up to the maximum full scale value.

Once phase 1 is finished, the message "Loaded weight - Waiting for stability" is displayed. Once the calibration value has been acquired, the zero acquisition phase begins and the display shows "Phase 2 - Unload weight" and the operator must proceed by removing the calibration weight from the pan.

The message "Unloaded Weight - Waiting for data computing" appears on the display and if all the phases are respected and the weights are correct, the calibration is accepted, the display shows the message "Instrument Calibrated" for a few seconds and then returns to the screen weighing standards.

If, on the contrary, errors have occurred, the display shows the message "Calibration Error" for a few seconds, then returns to the standard weighing screen and the previous calibration value will be kept.

**NOTE: The calibration process is immediately interrupted if there are more than twenty grams on the weighing pan at the time of calibration.**

## 10 QUICK RESTORE OF DEFAULT PARAMETERS

If you have mistakenly changed the SETUP parameters with an invalid configuration, you can restore the default parameters (factory settings) through a simple method.

- Remove the power supply connector of the scale located on the back of the same.
- Press and hold the first multifunction button on the left.
- Reinsert the power connector
- Wait for the word "SET FACTORY PARAMETER" to appear on the display.
- Now release the previously pressed key.
- In this phase the scale automatically restores the factory parameters

Note:

When a configuration has been modified and is correct, it is advisable to immediately use the Set Default

function in the SETUP menu (see 9.3.10) in order not to lose (if the instrument is turned off) your own personalized configuration.

## 11 SERIAL OUTPUT

EU-C Gold serial output is compatible with EIA RS232 standard.

Every answer from balance is sent every 100 ms.

### 11.1 SERIAL OUTPUT CONNECTION

BALANCE	DB9 CONNECTOR (9 pole)	DB25 CONNECTOR (25 pole)
RS 232 Out	Pin 2	Pin 3
RS 232 In	Pin 3	Pin 2
GROUND	Pin 5	Pin 7

### 11.2 DATA TRANSMISSION FORMAT

The format of the data transmitted by the balance is different as a function of the selected protocol by the menu option 9.2.1.3.

EURO and CRYSTAL protocols are suitable for computer interfacing; PRINTER protocol may be used to drive external printers.

#### AVAILABLE PROTOCOL TYPES

##### 11.2.1 CRYSTAL protocol

The answer to any type of command is formatted with a data field, a unit field and a status field as follows:

a measure field

a unit field

a status field

D0	D1	D2	D3	D4	D5	D6	D7	D8	D9	B	U	U	U	B	F1	F2	CR	LF
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18	19

The measure field (1..10) is the reading on the display.

It is right justified with decimal point and measure sign.

Unit field shows the current unit weight

Status field show the current balance status

B is the blank character.

Field N°18 and 19 are:

**CR** carriage returns ASCII code.

**LF** line feed ASCII code.

#### Status information of the F1 character

<b>D</b>	Transmitted data is valid
<b>O</b>	Balance in over range
<b>U</b>	Balance in under range
<b>T</b>	Tare
<b>C</b>	Calibration process in progress
<b>Z</b>	Zero acquisition
<b>I</b>	Init test

**Status information of the F2 character**

<b>S</b>	Measure is stable
<b>I</b>	Measure is not stable
<b>E</b>	Error
<b>A</b>	Automatic calibration (when F1 = C)
<b>L</b>	Load weight(manual calibration) (when F1 = C)
<b>U</b>	UnLoad weight (manual calibration) (when F1 = C)
<b>D</b>	Calibration done (when F1 = C)
<b>B</b>	Calibration busy(already started)
<b>O</b>	Calibration OFF
<b>P</b>	Percentage

**11.2.2 EURO protocol**

The answer to any type of command is formatted as follow:

D0	D1	D2	D3	D4	D5	D6	D7	D8	CR	LF
1	2	3	4	5	6	7	8	9	10	11

The measure field (1..9) is the display reading.

It is right justified with decimal point and the measure sign.

Field N°10 and 11 are:

**CR** carriage returns ASCII code.

**LF** line feed ASCII code.

**11.2.3 PRINTER protocol**

Is activated by the "PRINT" pushbutton. The output string uses 10 characters for the measure including sign and decimal point, the value is right justified and filled with spaces from the left. The format of the output is:

D0	D1	D2	D3	D4	D5	D6	D7	D8	D9	B	U	U	U	CR	LF
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16

The measure field (1..10) is the display reading.

It is right justified and has the decimal point and the measure sign.

Unit field shows the current unit weight

B is the blank character.

Field N°15 and 16 are:

**CR** carriage returns ASCII code.

**LF** line feed ASCII code.

**Note**

Data will be printed if and only if the weight on pan is stable.

**11.3 SERIAL COMMANDS**

The balance responds to the following commands only if a CRYSTAL or EURO protocol is selected:

SERIAL REMOTE CONTROLS TABLE

COMMAND	FUNCTION
<b>B</b>	Send data on display
<b>C</b>	Start calibration
<b>I</b>	Start continuous transmission of data on display
<b>F</b>	Stop continuous transmission of data on display
<b>T</b>	Tare command

**Note 1:**

Every remote control command must end with <CR>.

Line-feed characters are ignored.

customer care

info@gibertini.com

Tel. (+39) 023541434

ISO 9001:2015

LAT N° 094

**GIBERTINI  
ELETTRONICA s.r.l.**

Via Bellini 37  
20026 Novate Mil.se  
Milano (Italy)

Tel. (+39) 02 3541434  
Fax (+39) 02 3541438  
www.gibertini.com  
sales@gibertini.com